

FIBER OPTIC ILLUMINATOR INSTRUCTION

ATTENTION CHANGE THE LIGHT BULB

Press the button of the top, and slide the top cover to the back.

STOP !

You have to **UNPLUG THE POWER CABLE** before open the box! and let the **LIGHT BULB COOL DOWN**



Remove the top of the cover and open the box.

Take the light bulb out off the housing.



Change the new light bulb and put it back.

Put the top cover back.



CHANGE THE SOCKET

Use the screw driver to loose the two screws showing in the picture, then you can take the socket out, and change the new one



HOW CAN ATTACHED THE ILLUMINATOR AND THE FIBER OPTIC CABLE

Push the adapter cap of the illuminator nose piece, and plug the cable adapter in.



When use the double pipe light guide, after you plug in the cable adapter, you need to use the screw driver to tight the two screws on the side of the nose piece of the illuminator



Warranty information

The light source for one (1) year warranty, the light bulb is not warranted.

Howard Electronic Instruments, Inc. Address: 974 SE Pioneer Road. El Dorado, Kansas 67042

Tel: 1-800-394-1984 Fax: 316-321-2803 Website: www.heiscope.com Email: sales@heiscope.com