**General instruction for STDO21 electrode**

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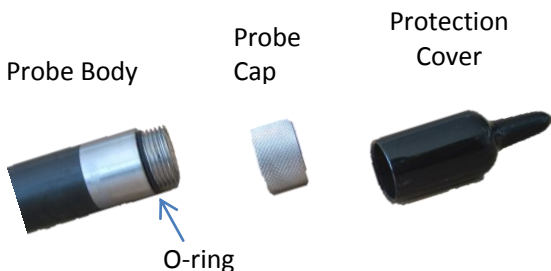
**Specifications:**

Connection:	Min-Din
Shaft length:	125mm
Shaft diameter:	13mm
Shaft material:	ABS
Temperature range:	0-60°C
Measurement range:	0.02~20 mg/L
Accuracy:	0.02mg/L

This STDO21 is a luminescence DO probe which need very low maintenance. Please perform calibration before first usage; then the probe does not calibration in a few months. We recommend customer to replace the **probe cap** about **every 12 months** or when STDO21 performance is not good any more.

When customer replaces a new probe cap, you **MUST** input the new probe cap parameter into the meter immediately, then do calibration before any usage.


**Replace Probe Cap Instructions**

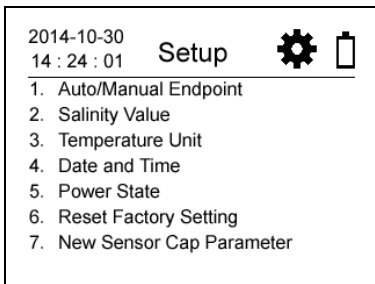30253566	DO probe cap, STDO21, 1 unit
----------	------------------------------


For optimum performance, use the following operating procedures:

1. Take off the old probe cap: One hand holds probe body, the other hand slightly turns probe cap counterclockwise until it's completely off from probe body.
2. Check whether O-ring is damaged: Please remove the old O-ring if there is any damage and replace a new O-ring.
3. Please make sure the new O-ring is clean. Please use clean cloth to wipe out any dusts or stains.
4. Put on new probe cap: one hand holds the probe body, the other hand gently turn the probe cap clockwise until the cap is tightly secured.
5. Connect STDO21 to ST400D meter.
6. Input the probe cap parameter

To enter the parameter input setting, you need to press READ + CAL + Setup button together when "7. New Probe Cap Parameter" is selected.

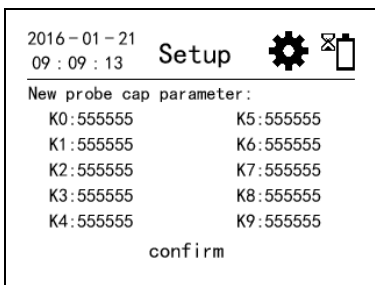Then meter goes to parameter input setting like following, input the parameter of this probe cap:

7. Calibration: Do 1 point or 2 points calibration before any usage.

## Storage

Rinse or clean the electrode after each use. Then keep it in a moisture and clean environment. Make sure the sponge in the protection cover is moisture.

## Precautions and Limitations

1. Do not expose the shaft to organic solvents, when cleaning or when taking measurements.

2. STDO21 has a built-in temperature probe; do not use the probe outside the recommended temperature range.
3. STDO21 can only work with ST400D meter.

Modelo	Nº de ítem	Descripción
STDO21	30219117	Sonda Óptica DO STDO21, 1.0 m cable
STDO21	30276108	Sonda Óptica DO STDO21, 5.0 m cable

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Ohaus Corporation, all rights reserved/ todos